

INSTITUTE FOR SUSTAINABLE TECHNOLOGIES NATIONAL RESEARCH INSTITUTE – RADOM

PW-004 Multi-Year Programme

DEVELOPMENT OF INNOVATIVE SYSTEMS OF MANUFACTURING AND MAINTENANCE 2004–2008




T-17 RECIPROCATING TESTING MACHINE WITH PIN-ON-PLATE TRIBOSYSTEM FOR TRIBOTESTING OF BIOMATERIALS




MAIN CHARACTERISTICS

T-17 Testing Machine is intended for the evaluation of combinations of sliding materials (e.g. metal-polymer couples) to be used as bearing surfaces of total-joint replacement prostheses for humans. The research is carried out on reciprocating motion.

Experiments can be carried out according to **ASTM F 732**.

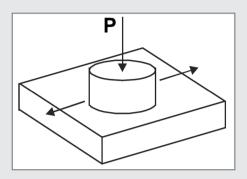

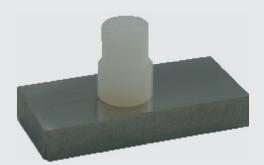

The tribosystem consists of the stationary test pin and the plate moving at the desired oscillatory frequency and amplitude. The test pin is pressed against the plate at the defined load P. The lubricant may be optionally delivered to the contact zone in a circulating lubrication system equipped with a peristaltic pump, heating device and filter.

T-17 Testing Machine is equipped with a control-measuring system that consists of the following:

- A set of measuring transducers,
- Controller,
- Digital measuring amplifier,
- PC and special software for measurements and data acquisition,
- Peristaltic pump (option), and
- Lubricant heating and filtering system (option).

During the tests, the following quantities are measured:

- Friction force.
- Total linear wear of test specimens,
- Ambient temperature,
- Lubricant temperature (option), and
- Time and the number of cycles (sliding distance).

Before the run, the desired frequency and amplitude of reciprocating movement are set. Then, the measured values are displayed on the monitor screen and saved on the computer disk. The motor of the tribotester is automatically stopped when the preset sliding distance (number of cycles) is reached. After test completion, one can print a report presenting the curves of the changes in the particular quantities versus time.

TECHNICAL SPECIFICATIONS

Type of movementContact geometry

Nominal pin diameter

Nominal plate dimensions

Amplitude

Frequency

Normal load

Lubricant temperature

• Tribotester dimensions (W x H x D)

Tribotester weight

Power supply

Max. power consumption

sliding - reciprocating

conformal

9 mm

 $36 \times 17 \text{ mm}$ (rectangular prism) or $\Phi 36 \text{ mm}$ (disk)

0.5-12.5 mm

0.5-9.5 Hz

up to 245 N

37°C (stabilised)

600 x 430 x 200 mm

30 kc

230 V / 50 Hz (optionally 110 V / 60 Hz)

0.3 kW

